



Evolve 128 Camera Update v9.13

(Released December 16, 2011)

Release notes:

The camera update application will update the following features for your Evolve 128 camera:

- Trigger First Mode with updated Overlap functionality provides higher frame rates.
- Strobe Mode Triggering with updated Overlap functionality provides higher frame rates.
NOTE: In overlapped Strobe mode, the camera will ignore Shutter-Open Delays
- Improved Dual-Camera triggering performance.

To use the higher frame rates provided by the updated triggering behavior, ensure that the camera is in “Pre-Sequence-Clear Mode” when acquiring images.

Usage instructions:

1. Download the camera update application.
2. Connect and turn-on only the Evolve 128 you intend to update.
3. Run the camera update application. *Do not power off or disconnect the Evolve 128 while running the camera update application!*
4. Accept the prompt to update – the application will determine your Evolve 128 camera version and apply the update accordingly.
5. After the update has completed, be sure to power-cycle the camera to ensure that all updates are applied.

If you have any questions or require further assistance, please call Photometrics Customer Service and Technical Support at +1 800.874.9789 or email support@photometrics.com.

HIGH PERFORMANCE EMCCD & CCD CAMERAS FOR LIFE SCIENCES