

Windows® Installation Notes for PVCAM® 2.6.9.3

PVCAM version 2.6.9.3 has been modified to address an issue with installation on computers running Windows XP. In addition, the ICL library has been enhanced with new functionality.

The installer will prompt you to update the firmware located on the PCI interface card, if required. You **must** upgrade to the latest version of the PCI firmware for this PVCAM version and the device driver to work correctly.

Installing PVCAM 2.6.9.3 on a clean system

To install PVCAM 2.6.9.3 on a system that has not had a previous installation of PVCAM, follow these instructions:

1. Install the PCI card!
2. Run the PVCAM 2.6.9.3 installation program (PVCAMSetup.exe) from the CD-ROM or hard drive. Choose to exit the installation without rebooting.
3. Restart the computer. You can run RSConfig if you want to change the default name of the camera, but this is not necessary.
4. Run PVCAMtest (found under the Start menu at Start>Programs>Roper Scientific>PVCAMtest) to verify that everything is working correctly.
5. (Optional) Run RSImageSetup.exe on the CD-ROM or hard drive to install the RS Image™ application.

Upgrading PVCAM when an older version was previously installed

To check if an earlier version of PVCAM has been installed on your computer, do a search of your hard disk(s) for the file PVCAM32.dll.

1. Run PVCAM 2.6.9.3 and select the Remove option to uninstall the older version.
2. Run the installer and install the new version following the instructions presented.
3. Run PVCAMtest to verify that everything is working correctly.