

## Free Camera Trial Pre-Arrival Guide

Thank you for your interest in a free trial of one of our Teledyne Photometrics products. We hope that our free trial process provides you with an appropriate time for evaluation. Please use this guide to ensure that you're ready for when the camera arrives.

### Mounting the camera:

You will need a 1-inch C-mount adapter at 1x magnification to mount the camera onto the microscope. Please ensure that the adapter has no additional lenses or magnification as our products will perform best without these.

### Using a suitable PC:

Our products are compatible with all Windows PCs/laptops running Windows 7, 8 or 10 – we recommend the 64-bit versions. You will need a minimum of 8GB of RAM, however we would recommend more (16GB+) for optimum performance.

All of our Teledyne Photometrics products connect through USB 3.0.

### Software and driver downloads:

We would recommend downloading our PVCAM camera drivers and Ocular imaging software prior to your loan. These can be downloaded from our website's [download page](#).

If you have any further questions, you can contact us through the live chat on our website, or your local sales representative.