

QIMAGING QUICK INSTALLATION GUIDE



Install the Camera-Interface Software

1. Power on the computer and insert the QImaging USB stick into your USB drive.
2. Install camera drivers by choosing "QI_setup.exe".
3. Follow the installation instructions on screen.
4. Install the USB3 PCIe card supplied with the camera.
5. Restart your computer and connect the camera to the USB port on the installed card.

Retiga™ 3000 and 6000



1. Using a USB "A to B" cable, plug the USB "B" end of the cable into the camera's USB "B" socket.
2. Plug the USB "A" end of the USB "A to B" cable into a USB port of your computer.

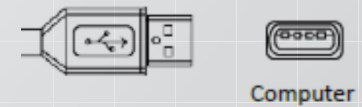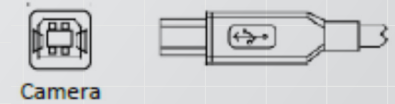

More detailed installation instruction can be found in the camera instruction manual on the USB stick.

Contact QImaging Customer Service for additional information: 1-800-874-9789 or www.qimaging.com



www.qimaging.com

©2015 QImaging. All rights reserved.

58-557-003 Rev A3