

QIMAGING QUICK INSTALLATION GUIDE


Install the Camera-Interface Software

1. Power on the computer and insert the QImaging USB stick into your USB drive.
2. Install camera drivers by choosing "QI_setup.exe".
3. Follow the installation instructions on screen.
4. Restart your computer and connect the camera to the PC.
5. If you purchased the QCapture Pro software, please follow the installation instructions on the QImaging USB in the Document folder.

Rolera™ Bolt



1. Using a USB "A to B" cable, plug the USB "B" end of the cable into the camera's USB "B" socket.
2. Plug the USB "A" end of the USB "A to B" cable into a USB port of your computer.

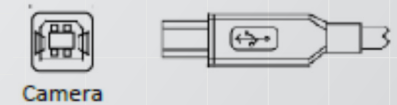

More detailed installation instruction can be found in the camera instruction manual on the USB stick.

Contact QImaging Customer Service for additional information: 1-800-874-9789 or www.qimaging.com



www.qimaging.com

©2015 QImaging. All rights reserved.

03-0023 Rev A2