

QIMAGING QUICK INSTALLATION GUIDE

Install the Camera-Interface Software

1. Install the high-speed PCIe card into the computer.
2. Power on the computer and insert the QImaging USB stick into your USB port.
3. Install PVCAM®.
4. Follow the installation instructions on screen.
5. Restart your computer and connect the camera to the host computer.

Install Ocular (If Available)


1. Insert the Ocular USB key into another USB port on the host computer. *This does not need to be a Super Speed USB3 port for proper operation. Wait for the Windows new hardware installation wizard.*
2. Run the SetupOcular installer file found on the QImaging USB stick. Note: the actual name will be "SetupOcular" followed by the Ocular version number.
3. Open the Ocular software from the Windows Start Menu.

optiMOS™



optiMOS™

1. Plug one end of the data cable into a high speed PCIe port of your computer. Plug the other end of the cable into the camera's data socket.
2. Connect the power cord to the power supply, plug the power cord into a wall outlet, and connect the power supply to the camera.
3. Power on your computer. Then power on the camera by pushing the power button.
NOTE: "WAIT" LED should be off before running the camera. It takes approximately 30 seconds for the camera to initialize.

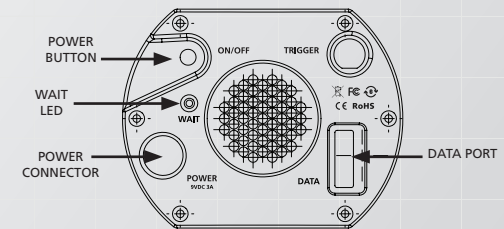

More detailed installation instruction can be found in the camera instruction manual on the QImaging USB stick.

Contact QImaging Customer Service for additional information: 1-800-874-9789 or www.qimaging.com



www.qimaging.com

©2017 QImaging. All rights reserved.

58-607-001 Rev A02